

Scanning Magnetic Microscope

The Tristan model 770 Scanning Magnetic Microscope is designed to measure magnetic fields above a planar surface with unparalleled spatial resolutions.

Using a liquid nitrogen SQUID sensor, it can be used to image diverse objects such as:

- shorts on ground planes
- traces on a circuit board or multi-chip module
- current distributions
- magnetic inks used in currency



Magnetic image of dollar bill section

- insulators, ferrous and non-ferrous metals to detect cracks, voids and corrosion
- nanoparticle distributions
- flux-motion in HTS materials


Besides measuring magnetic fields, the 770 can also be configured to detect:

- induced magnetization
- magnetic susceptibility
- eddy currents
- magnetic hysteresis
- Barkhausen effect
- micropaleontology
- magnetobiologic activity




Model 770 Scanning SQUID Microscope with dc and ac biasing magnets

- The 770 is a fully featured measurement system that allows the user to extract a magnetic image of the object being measured over the entire dc – 10 kHz frequency range.
- Whether the samples are circuit boards, multi-chip modules, steel or aluminum plates, composites or even plastics, the 770 can measure surface and even deeply embedded sources with a spatial resolution down to 50 μm .
- The use of a High Temperature Superconducting YBCO step-edge dc SQUID sensor gives it unparalleled sensitivity with the ability to measure fields smaller than 20 $\text{pT}/\sqrt{\text{Hz}}$. Tristan's HTS sensors can also operate in applied magnetic fields exceeding 1000 oersteds.
- The model 770 allows computer controlled scans of objects over a large (15 x 15 cm) area with 25 μm stepping capability with sub-micron stepping available. The user has the ability to preprogram the scan coordinates.
- The 770 requires minimal setup. Automated setup and computer control makes measurements rapid and repeatable. System software provides the ability to control the critical system components, acquire data from the SQUID sensor, and analyze the data to determine the magnetic properties of the sample being measured. The use of open architecture software allows the user to modify and customize nearly all aspects of operating including image processing.

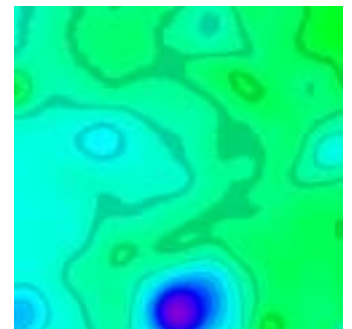

scan of concrete showing magnetic impurity

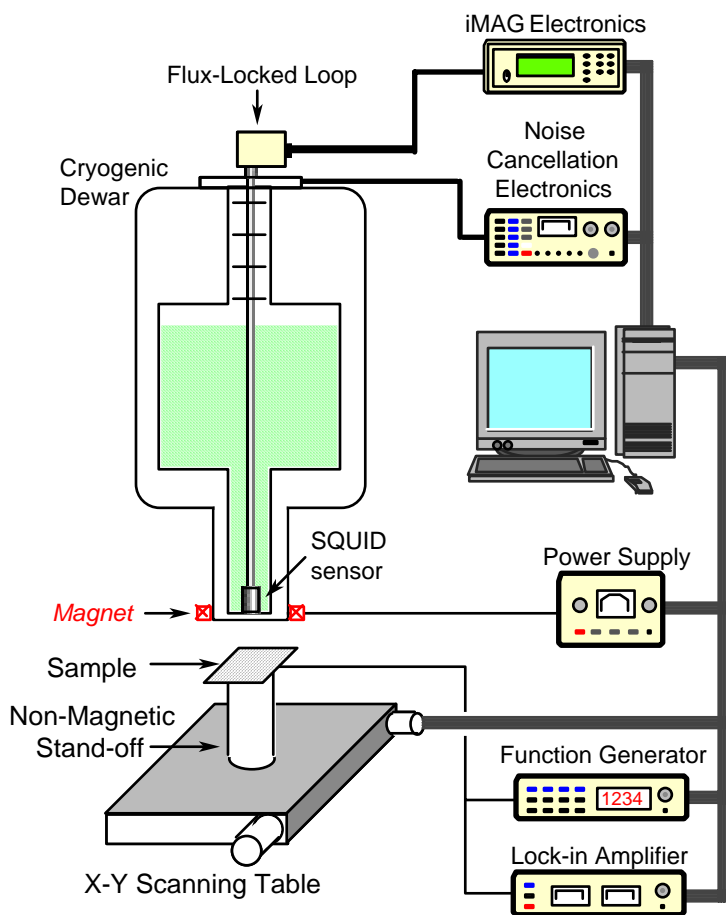

Components of the Model 770 system including optional accessories

The Standard model 770 is configured to detect electric currents and to measure remnant magnetic fields. It includes a Single-Channel Scanning SQUID Magnetometer Probe, iMAG[®] SQUID Electronics, Cryogenic dewar, Room Temperature Scanning Stage, Computer Control and Data Acquisition System, and Imaging Software. The model 770 can be supplied with additional capabilities to extend its measurement capabilities.

OPTIONS AND ACCESSORIES

SCAN AREA: Larger scan areas (e.g., 30 cm x 30 cm) and higher resolution stepping (25 μm standard) are available upon request.

Substitution OF HTM-1 SQUID Sensor: For measurements where sensitivities significantly below 20 pT/ $\sqrt{\text{Hz}}$ are needed and ultimate spatial resolution is not as important, Tristan can substitute the model HTM-1 sensor with a significantly larger (1 mm x 1 mm) detection coil with a sensitivity better than 3 pT/ $\sqrt{\text{Hz}}$.

Fixed Field Capability: This option consists of a fixed field that generates a vertical (B_z) field on the sample. This allows magnetic susceptibility measurements on insulators, conductors and ferrous materials to be performed.

ac Field Capability: This option allows a small vertical ac magnetic field to be imposed on the sample. This capability is of interest when eddy current measurements are desired. This option can be used simultaneously with the dc Field Option for added flexibility in magnetic characterization.

Horizontal Field (B_x and B_y) Sheet Inducer: A horizontal field sheet inducer, which can apply an ac magnetic field parallel to the test surface, to induce a large extended eddy current in a desired orientation, can be used to image cracks or material loss deep in conductive (e.g., aluminum) structures.

ac Field Compensation Electronics: When imaging conductive materials, if an ac signal is directly coupled into the system, the resultant signal (from the field coils) may be much larger than the signal from the sample. To minimize this, Tristan can supply an ac Compensation system to null the ac signal in the detection coil and extract the induced signal in the object being measured.

SPECIFICATIONS

SENSOR: High temperature superconducting quantum interference device (SQUID) operating at 77 K

SPATIAL RESOLUTION: Better than 100 μm

SENSITIVITY: 2×10^{-11} tesla/ $\sqrt{\text{Hz}}$ (20 pT/ $\sqrt{\text{Hz}}$)

OPERATING BANDWIDTH: dc - 10 kHz. Measurements can be made at any frequency. Bandwidths above 20 kHz are available.

CRYOGENIC COOLING: To avoid low frequency noise below 200 Hz, the system uses liquid nitrogen to cool the sensor.


CRYOGENIC HOLD TIME: Time between refills of liquid nitrogen is typically 3 days

SAMPLE SCANNING RANGE: 15 cm x 15 cm in x-y directions

SCAN STEP SIZE: Adjustable with minimum step size of 25 μm .

SAMPLE PREPARATION: None required. Samples are measured at room temperature

POWER REQUIREMENTS: 100, 115 or 220 VAC, 50 or 60 Hz

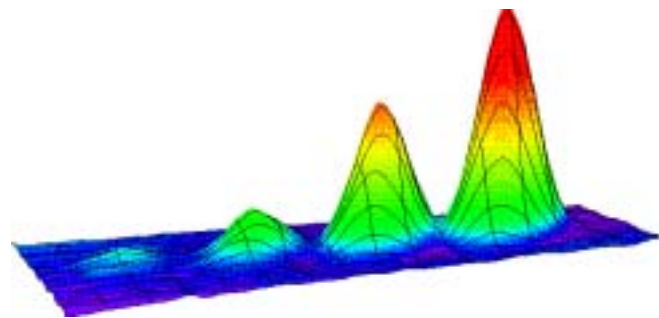

scan of holes in a steel plate



6350 Nancy Ridge Drive, Suite 102
San Diego, CA 92121
(858) 550-2700 [fax] 550-2799
<http://www.tristantech.com>



Represent by
NIKI GLASS CO., LTD. TOKYO

9-7 Mita 3-Chome Minato-ku Tokyo 108-0073 Japan
TEL : 81-3-3456-4700 FAX : 81-3-3456-3423